

QUICK GUIDE SMART

1. First start-up







During the first boot, the screen (*Figure 1*) will display proposing to save the "Download" folder (among others with instruction manual) on the USB flashdrive. Insert the USB flashdrive and wait a second to detect the hardware, then press . By pressing  you quit downloading the folder, the window will appear again during the next start. You can tick „Don't show again” so that the window will not be displayed when you start.

Figure 1 – Downloading the files




You can always download the "Download" folder in the information panel. Press , enter to info panel , press icon .

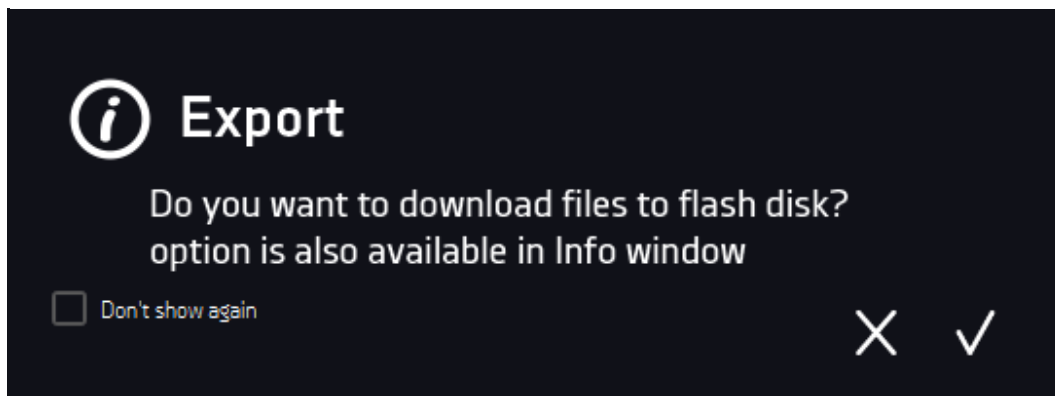
QUICK GUIDE SMART


2. Base window

The base window is used to read the current temperature in the chamber and to program and set the device parameters.

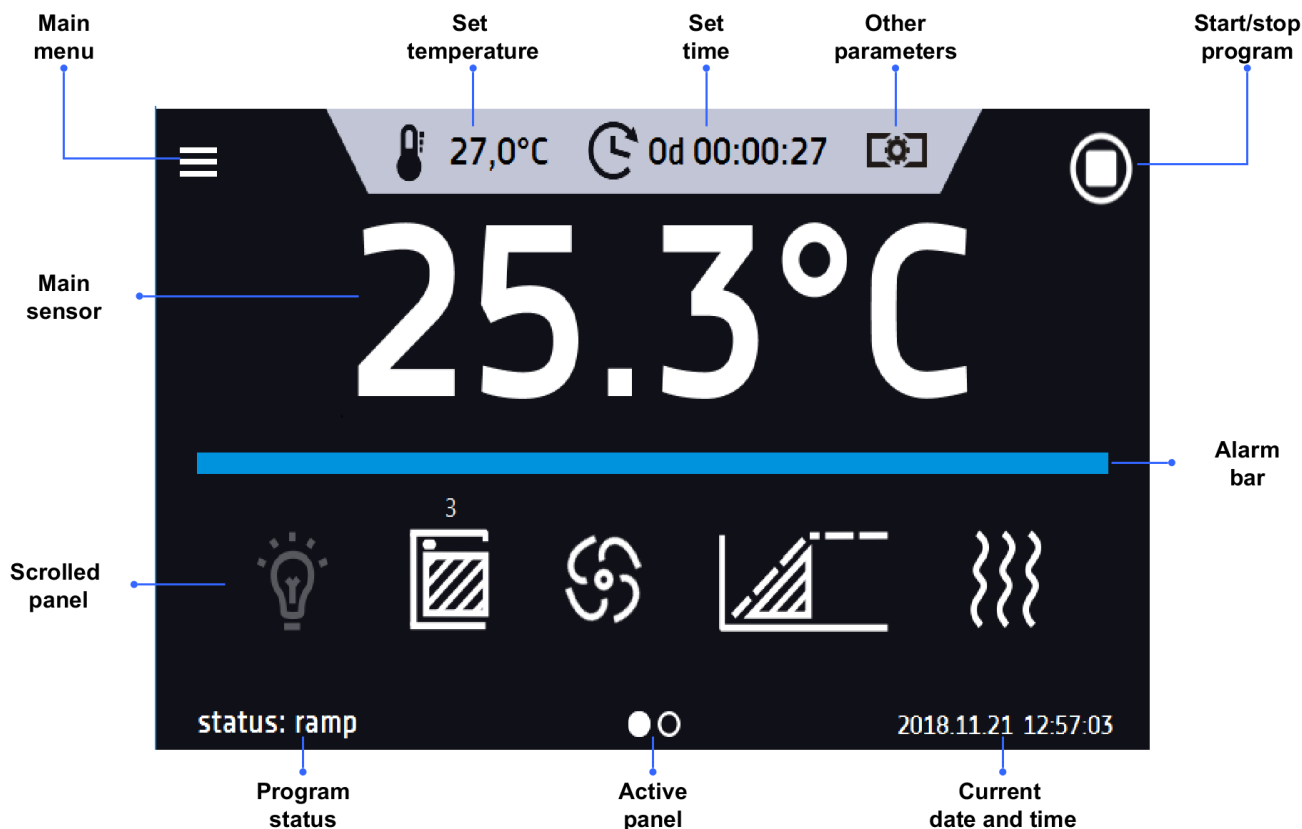


3. Quick Program





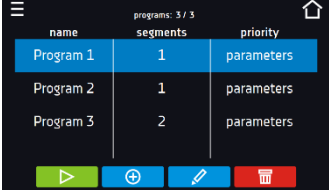

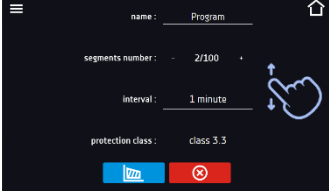


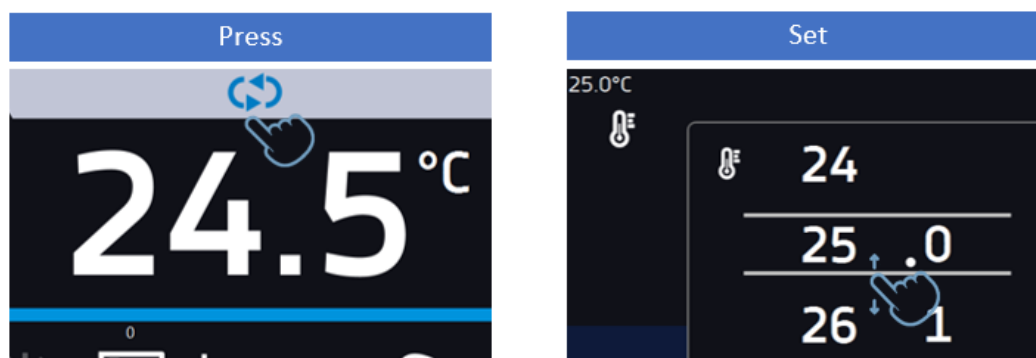Quick Program allows you to quickly start the program from the home screen position without having to enter and create a menu. After switching on the device, a symbol  will be available on the main screen.



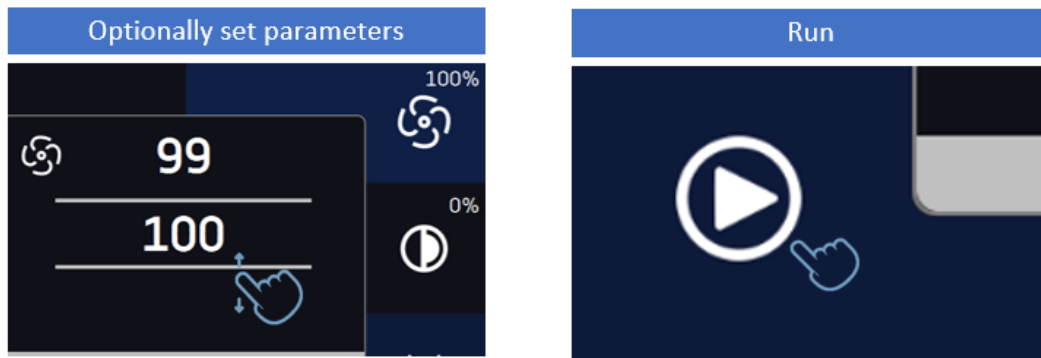

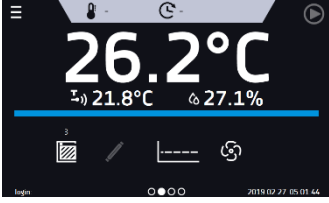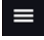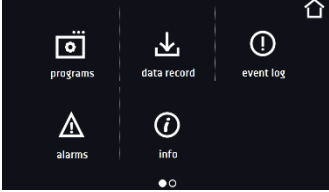


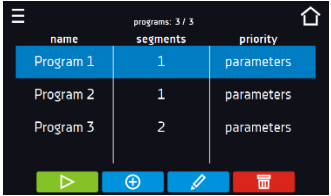


QUICK GUIDE SMART






4. CREATING AND STARTING UP THE PROGRAM

	<p>Press  in the base window to enter into menu.</p>
	<p>Press programs .</p>
	<p>In Programs panel press .</p>
	<p>Set the other program parameters (with more parameters you can scroll the panel up and down). Go to program segments by pressing .</p>
	<p>Using the side keyboard, set the parameters of the Temperature, Time (the number of parameters depends on the device and additional options). If you want the device to reach the set temperature as quickly as possible, allow the ramp time to 0.</p>

QUICK GUIDE SMART

	<p>Navigating between the program parameters, segments and summary is done by swiping up the segment with the finger or touch the icon  .</p>
	<p>If you have set parameters for all segments, move to the last window - Summary. In this window, all segments are visible along with the entered settings. If everything is correct, move the window down and press .</p>
	<p>The saved program will be automatically displayed on the list. To start the program press .</p> <p>To stop a running program press .</p>

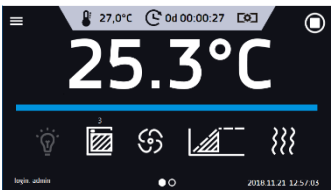5. QUICK START / STOP THE PROGRAM

	<p>In the main menu press:</p> <ul style="list-style-type: none">  - to start the program  - to stop the program
---	--